


DR Software Update Instructions

- 1) Choose the Software Download link for your instrument.
Example: If you have a DR2800 Spectrophotometer, click '[DR2800 Software Download](#)'

Note: *The download can take up to 1 minute.*

- 2) Click **Open**, when the following prompt is displayed: 'Do you want to **open** or save this file?'
- 3) Select all files ending with **.gz**

Note: *The number of .gz files varies between different instruments.*

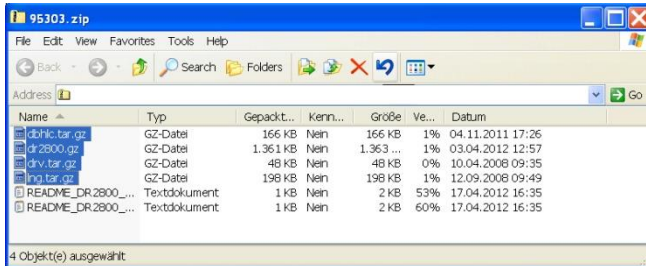

- 4) Copy all files with ending **.gz** to a USB flash drive.
Note: *Please do not unzip the single .gz files.*
- 5) Select '**System Checks**' in the main menu of your instrument. Then select '**Instrument Update**'.
- 6) Plug the USB flash drive into the USB port of your instrument.
- 7) Press **OK**. The software will be updated.
Note: *Please wait until the update is complete. This can take up to 10 minutes.*
- 8) When the instrument software has been updated, a prompt to restart the instrument is displayed. Switch off your instrument. Wait at least three seconds and switch on again.

Your spectrophotometer is now updated.